




IF YOU WANT TO:	THEN DO THIS:	DESCRIPTION
<p>Turn on/off power</p>	<p>Press the front, left I/O button.</p> <p>Press PAUSE to view the current settings and operate the oven.</p>	<p>Controls line power to the unit.</p> 
<p>Adjust the oven temperature</p>	<p>Press and hold the UP or DOWN button until the set point blinks to access the set point.</p> <p>Press the UP or DOWN button to set the temperature (Centigrade).</p>	<p>These controls are associated with the heating function.</p> 
<p>Adjust the rotisserie motor speed, pause the motor or jog the motor</p>	<p>Press and hold the UP or DOWN button until the set point blinks to access the set point.</p> <p>Press the UP or DOWN button to set the speed (RPM).</p> <p>Press PAUSE to pause the rotisserie.</p> <p>When the door is open, you can use the PAUSE button as the JOG button.</p> <p>Press PAUSE again to take the oven out of pause to program or run the oven.</p>	<p>These controls adjust the speed of the rotisserie motor.</p> 

[www.affymetrix.com](http://www.affymetrix.com) Please visit our web site for international distributor contact information.

# Quick Reference Card



## GeneChip® Hybridization Oven 645 Best Practices

- Never stack more than two ovens.



- Allow for adequate air flow, i.e., do not block rear vent.



THIS QUICK REFERENCE CARD WAS DEVELOPED TO AID IN THE PROPER USE OF THE AFFYMETRIX® GENECHIP® HYBRIDIZATION OVEN 645.

- Place ovens away from air conditioning units
- Ensure proper amount of free space around oven.



- Do not attempt to monitor the thermal performance of the oven using non-Affymetrix temperature probes. Affymetrix calibrates the temperature profile at the factory to ensure proper operation.



- Balance the carriers in the rotisserie by placing each carrier 180° from each other.



- Use the carrier retainer to secure the cartridge carriers.

